LCD Digital Microscope II

#44341

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LCD Digital Microscope II

Quick Start Guide- Model #44341

Thank you for purchasing the Celestron LCD Digital Microscope II. Your microscope is a precision optical instrument, made of the highest quality materials to ensure durability and long life. It is designed to give you a lifetime of enjoyment with a minimal amount of maintenance.

This Quick Start Guide will introduce you with the basic functions of the microscope. Please visit celestron.com to download the complete instruction manual. We recommend reading the manual to fully familiarize yourself with all the microscope's functions and operations.

This microscope provides magnification from 40x up to 400x (up to 1600x with digital zoom), ideally suited for examining specimen slides of all kinds. You will view specimens or objects on the LCD Monitor, instead of through eyepieces found on traditional microscopes. With the LCD Monitor, specimens or objects are easy to see, and can be viewed by multiple people at one time. You can also take snapshots or short videos with the built in digital camera.

ACCESSORIES INCLUDED WITH YOUR MICROSCOPE

- 2GB SD Card
- USB 2.0 Cable
- 5 Prepared Slides
- 4 plug, multi-country AC adapter
- Dust Cover
- Carrying Case

SPECIFICATIONS

Mechanical Stage	3.5" x 3.5" (88mm x 88mm)
Digital Camera	5 MP CMOS; 10x Magnification in lieu of an eyepiece; 3072 x 1728 pixel array
LCD Monitor	3.5" with 4x Digital Zoom – High definition digital TFT display Resolution – 320 x 240 pixels, Contrast ratio – 350, Color Support – 262K colors, Brightness/Luminance – 400cd/m ²
Focuser	Smooth, single motion
Memory	SD card slot, supports up to 32GB card - 2GB SD card included (1,250 snapshots at 5MP, 40 minutes of SD video)
USB Cable	1.5 m long, 2.0 USB
Filter Wheel	6 Position- Red, Blue, Green and 1, 3, 6mm DIA aperature
Nosepiece	3 Position with click stop
Illumination	Upper and lower, built-in adjustable LEDs
Condenser	N.A. 0.65
AC Adapter	Input Universal 100 to 240Volt 50/60HZ
Weight/ Dimensions	56.44 oz/1600 g
Batteries	4AA (User supplied- up to 3 hours of use)

MAGNIFICATION (POWER)

Use the following table to determine the magnification power combinations you can achieve with the LCD Digital Microscope II.

Objective Lens	4x	10x	40x
Magnification of image on screen	40x	100x	400x
Maximum with 4x digital zoom feature	160x	400x	1600x

WORKING DISTANCE

Objective Lens	4x	10x	40x
Working Distance (mm) - est.	35.3	7.8	0.7

SETTING UP YOUR MICROSCOPE

1. Remove the carrying case from the shipping box.
2. Unzip the case, and carefully remove the microscope and other parts and set them on on a flat surface.
3. Insert the cylindrical plug on the end of the AC adapter into the socket on the back of the microscope (Figure 2) and then connect the plug end of the AC adapter into a proper power outlet as shown in Figure 3.

Note: The AC Adapter's plug is interchangeable for use around the world. Switch easily among USA, UK and EU style plugs.

To change the plug, push down and hold the small button release (circled in Figure 4). Grasp the plug prongs with your free hand and rotate the plug head counterclockwise 90° to release. Reverse the procedure with your chosen plug head. When the new plug head is secured in place, you will hear a click.

Battery Operation – You can use your microscope without AC power if you choose. This gives you the freedom to operate the microscope outdoors or indoors anywhere you please. You'll need 4 AA batteries (user supplied). Open the battery door on the bottom of the microscope and insert the batteries according to the battery polarity shown in the battery compartment. After the batteries are installed, close the battery door. Battery life will typically be three hours.

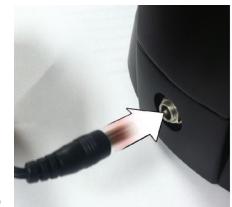
Fig 2

Fig 3

Fig 4

USING AN SD CARD

The LCD Digital Microscope II comes with a 2GB SD card, which you can use to capture images (snapshot or video). SD cards are inserted into the SD card slot in the LCD monitor (Figure 1).

Product design and specifications are subject to change without prior notification. This product is designed and intended for use by those 14 years of age and older.

© 2019 Celestron • All rights reserved • celestron.com/pages/technical-support • 2835 Columbia Street • Torrance, CA 90503 USA • 10-19 celestron.com/support/warranties

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