INFINIVIEW™

LCD DIGITAL MICROSCOPE

INSTRUCTION MANUAL

MODEL #44360, #44361

2. INTRODUCTION

Thank you for purchasing the Celestron InfiniView LCD Digital Microscope with a 5MP sensor and 3.5" color LCD screen. Your microscope is a precision optical instrument, made of the highest quality materials to ensure durability and long life. It is designed to give you a lifetime of pleasure with a minimal amount of maintenance.

Before attempting to use your microscope, please read through the instructions and reference the diagrams below. Familiarizing yourself with the many features of the Celestron InfiniView microscope will go a long way to making your experience enjoyable.

The InfiniView microscope provides low power viewing from 4x to 160x using the built-in software and LCD screen. For additional magnification up to 800x, connect the microscope to your PC. (Magnification depends on monitor size.) InfiniView is a low magnification power microscope, mainly suited for examining specimens like coins, stamps, rocks, electronics, insects, and other 3D objects. Specimens on slides can be viewed with top illumination, but this microscope is not ideal for this purpose. See some of our biological microscopes more suited to this type of viewing at celestron.com.

The Celestron InfiniView LCD Digital Microscope does not use an eyepiece like a traditional microscope. Instead, the camera sensor "sees" the specimen with 10x magnification, and the image is displayed on the high resolution color LCD screen. This makes viewing fun, cuts down on neck and eye strain and allows for easy sharing with others. Recording still images and video is a snap. Saving, transferring and sharing your image files is quick and easy with the included 2.0 USB cable and your home computer. Additionally, you can view on most TV screens with the AV/TV Cable.

Take some time to read the simple care, maintenance and troubleshooting tips provided for you in this manual to ensure that your InfiniView microscope provides you with years of quality performance, usage, and enjoyment.

IN THE BOX

InfiniView
LCD Digital
Microscope (1)

AC Adapter (1)

Lithium Battery (1)

USB cable (1)

TV Out cable (1)

QSG (1)


Cleaning Fabric (1)

Calibration Ruler (1)


PARTS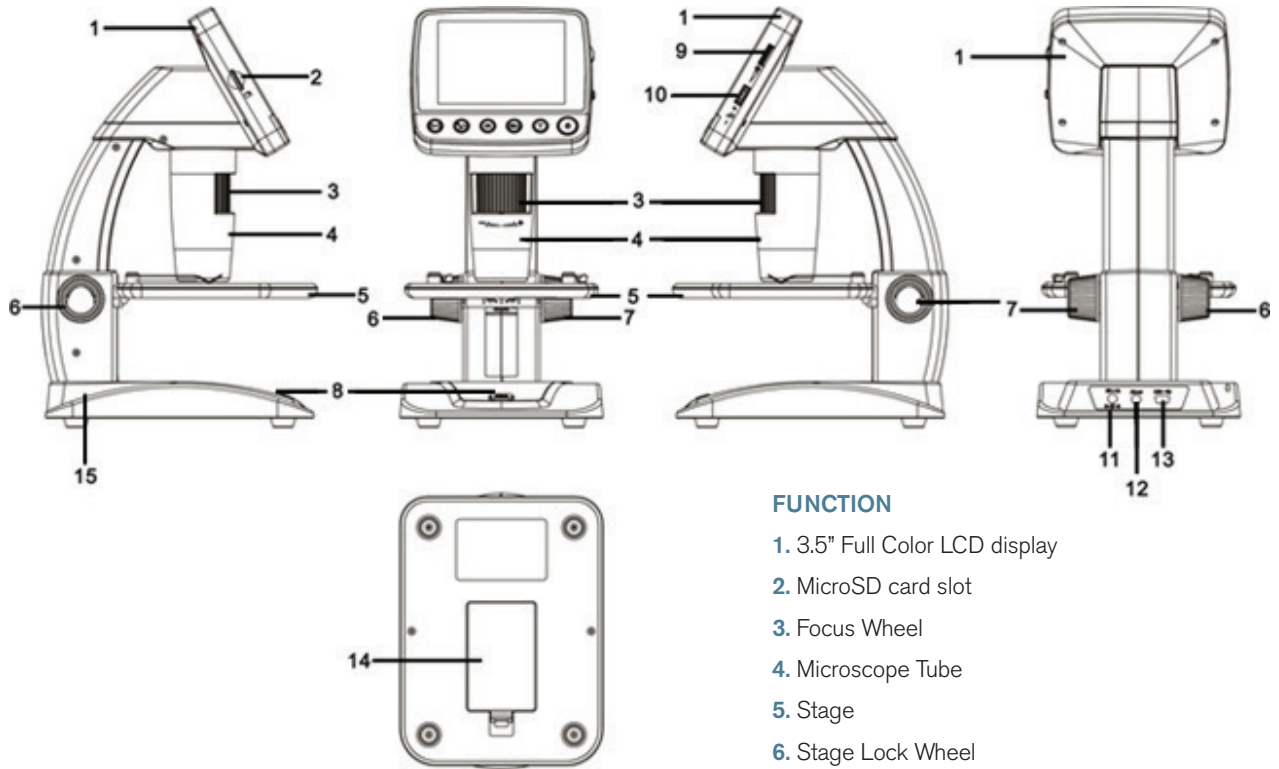

FUNCTION

1. 3.5" Full Color LCD display
2. MicroSD card slot
3. Focus Wheel
4. Microscope Tube
5. Stage
6. Stage Lock Wheel
7. Stage Height Adjustment Wheel
8. Capture Button
9. LED Illuminator Adjustment Wheel
10. Digital Zoom Switch (+/-)
11. 5v DC In port
12. TV Out port
13. USB port
14. Battery Compartment
15. Battery Charge Indicator LED

LCD SCREEN DETAIL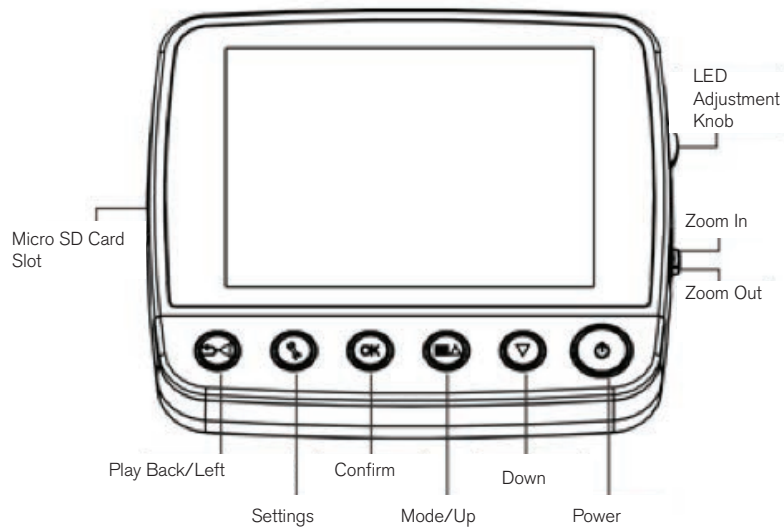

SPECIFICATIONS

SKU	44360 and 44361
LCD Monitor	3.5" (90mm) TFT Color LCD /4X digital zoom/320 x 240 pixels, 4:3 ratio
Sensor Size	5 MP (mega pixels)
Stage Adjustable Height	4.38" x 4.38" x .38" (111mm x 111mm x 9.5mm)
Lens	5 element IR cut , high quality glass/EFL(Equivalent Focal Length): 15.8mm /FOV: 13.0°
Magnification	4x to 160x using LCD screen (Up to 800x on computer monitor via USB connection)
Focuser Range	10mm to infinity
Image Distortion	.10%
Illuminator	Upper ring- 8 bright white adjustable LEDs
Shutter Trigger	On base – limits image shake
Still Image Resolution	Adjustable: 5MP, 3MP, 1.3MP, VGA (12MP with interpolation*)
Video Resolution	VGA, 30fps
Power	800 mAh rechargeable lithium battery
Memory	MicroSD card slot- Supports up to 32G (card NOT included)
Computer Interface	2.0 USB- Battery charging option through USB also
AC Adapter	44360: US Plug 44361: (multi country plug)- 100 to 240V- 50/60Hz
Software	On unit: Imaging software with zoom and timer features Downloadable from Celestron.com: Celestron Portable CapturePro (with measuring and note-taking features)
Operating Systems Supported	Windows: Windows XP/Vista/7.0/8.0 Mac: OSX 10.6/10.7
Software Languages	On unit: (9) English, German, French, Spanish, Italian, Portuguese, Japanese, Chinese Simple, Chinese traditional On Computer: (12) English, German, French, Spanish, Russian, Italian, Portuguese, Dutch, Polish, Japanese, Korean and Chinese
TV Output	AV cable (included)
Dimensions	149mm x 104mm x 227mm (5.80" x 4.10" x 8.90")
Weight	1.5 lb /24 oz. or 680 grams


*Interpolation refers to software adding pixels to an image to artificially make it larger. This can be beneficial when printing in larger formats.

3. SETUP

INSTALL BATTERY

Open the Battery Compartment and insert the Lithium battery, making sure to align the connection points correctly.

NOTE: Although the battery has a partial charge, if you are going to use the microscope with the battery only, be sure to charge the battery first. See below for more details on charging the battery.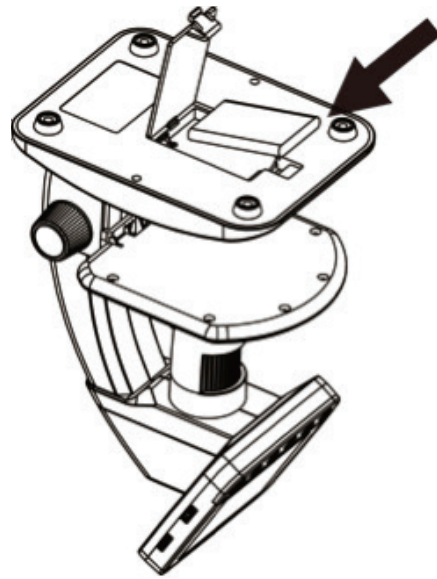

INSERT MICRO SD CARD

To capture images with your microscope while NOT connected to a computer, you need to insert a MicroSD card.

NOTE: The MicroSD card is NOT included. To capture images without a MicroSD card, see section 5.E in this manual for instructions on installing the computer software and using it to save images to your computer.

Locate the MicroSD card slot (2) on the side of the LCD housing.

Slide the card into the slot until it clicks into place and latches. Be sure not to force the card into the slot, as this may damage the card or the slot.

NOTE: The MicroSD card should be formatted before use. See section 6.2.G for details on formatting the card.


POWER ON



Press the POWER button to turn the unit on.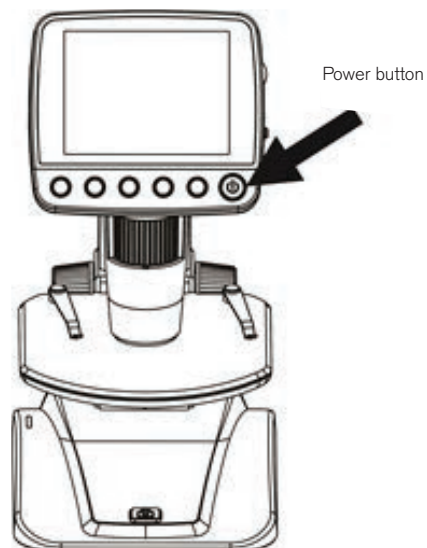


Press POWER button again to turn the unit off.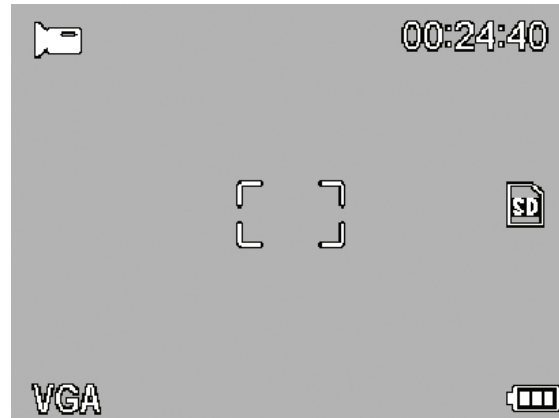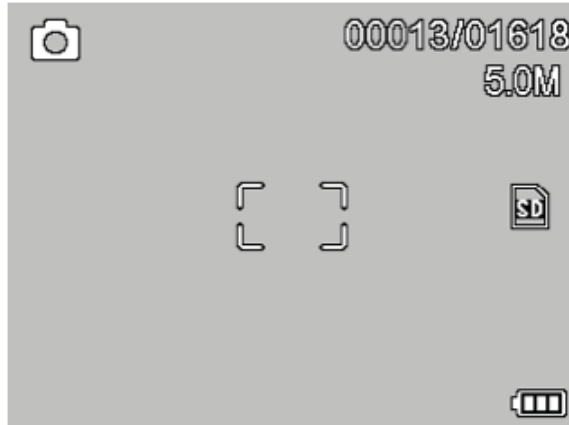

NOTE: The InfiniView has a Power Save feature. Refer to section 6.2.C for details.

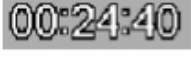DISPLAY ICONS

	Small Image Capture
	Video Capture
	Time Lapse Capture
	Number of Images Taken/Number of Images Remaining
	Video Capture Time Remaining (Hours:Minutes:Seconds)


	MicroSD Card Recognized
	Video Capture Setting
	Current Resolution Setting
	Battery Power Indicator

NOTE: The Display Icons can be turned off by pressing the Zoom Button (10) inward.

CHARGE THE BATTERY


- + Plug the small end of the power adapter into the DC In port (11) located on the back of the unit.
- + Plug the other end into a standard AC wall outlet.
- + The Battery Charge Indicator LED (15) will light up red to show the battery is charging.
- + When battery is fully charged, the Battery Charge Indicator LED will turn off and you can unplug the unit.
- + You can now use InfiniView without an external power source continuously for up to 2 hours.
- + You can monitor the battery power level on the LCD screen. Look for the Battery Power Indicator Icon described above.

NOTE: If you purchased your InfiniView outside of North America (item 44361), you will notice that the AC adapter has 3 interchangeable plug heads. Refer to section 9 to see how to change the plug heads.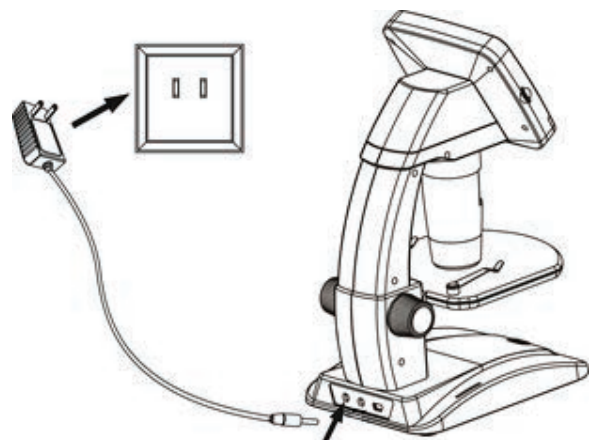


4. OBSERVING AND CAPTURING IMAGES: DIGITAL IMAGING

A. Make sure your InfiniView is powered ON.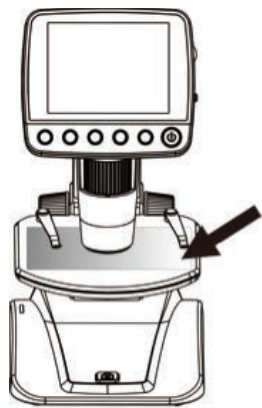


Put object onto the stage

B. Place your specimen on the stage (5). If your object is thin enough, secure it to the stage using the chrome clips.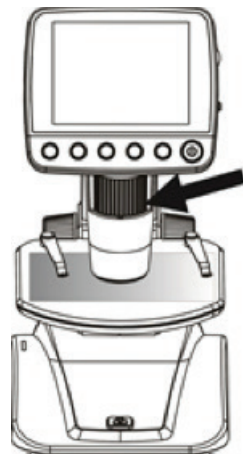

D. Rotate the Focus Wheel (3) to get a fine focus on the object.

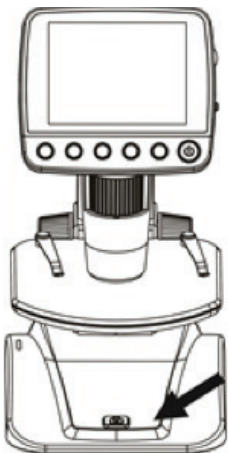
F. To record an image, press the Capture Button (8).


NOTE: Press the Playback/ Left button on the LCD screen housing to scroll between Still Image Capture mode and Video Capture mode.

C. Rotate the Stage Height Adjustment Wheel (7) to bring your specimen closer to the Microscope Tube (4).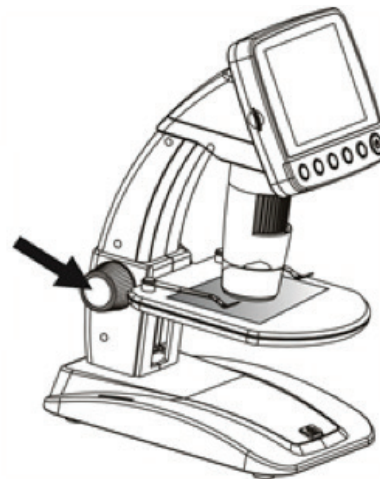

E. When the desired focus is achieved, lock the stage in position by tightening the Stage Lock Wheel (6).


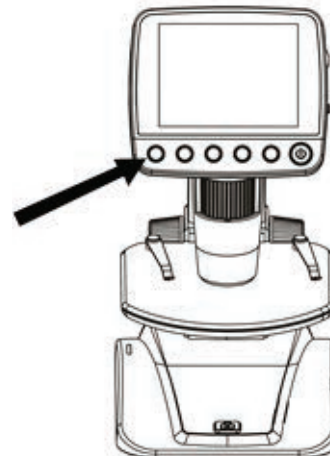
G. While in video mode, press the Capture Button to start recording video. Press the button again to stop.

5. SOFTWARE FEATURES

REVIEW SAVED IMAGE FILES:

To play back video or review a still image, simply press and hold the Playback/Left button on the LCD housing.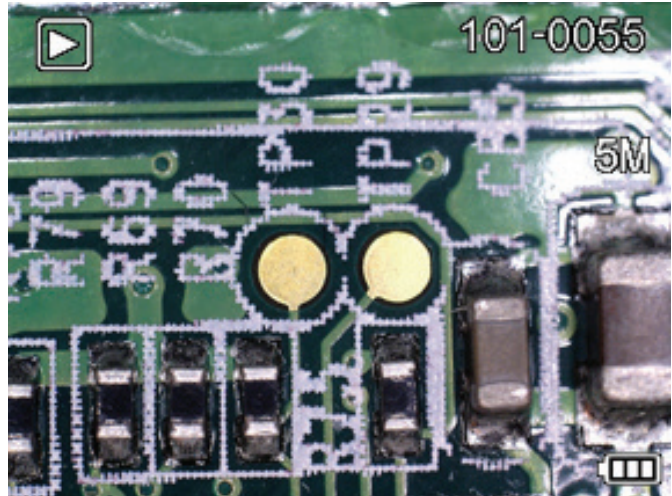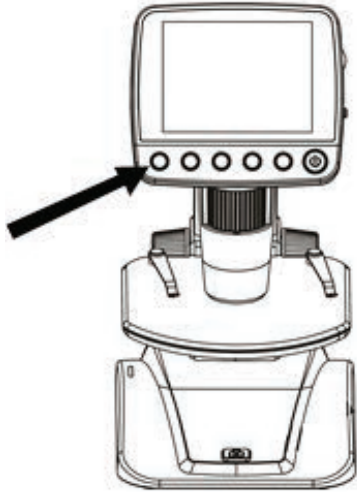

The latest image (or a thumbnail of the latest video) will appear on screen, with a directional arrow at the top of the screen.

To scroll through the images and thumbnails, press the UP  and/or DOWN  button.


To modify a particular image or video using the Playback feature, press the Settings button. 

The following screen will appear with 3 choices:

1. Delete
2. Protect
3. Slide Show


1. DELETE IMAGE FILES:

This choice deletes the selected image/video or ALL images.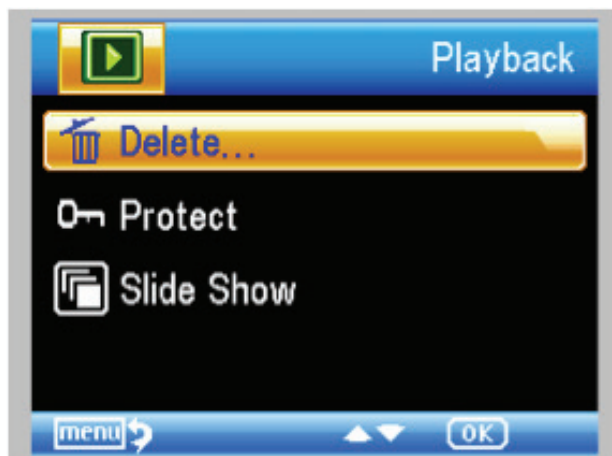


2. PROTECT IMAGE FILES:

This choice allows you to 'lock' and 'unlock' the selected image/video file or ALL images/video files on disk for protection.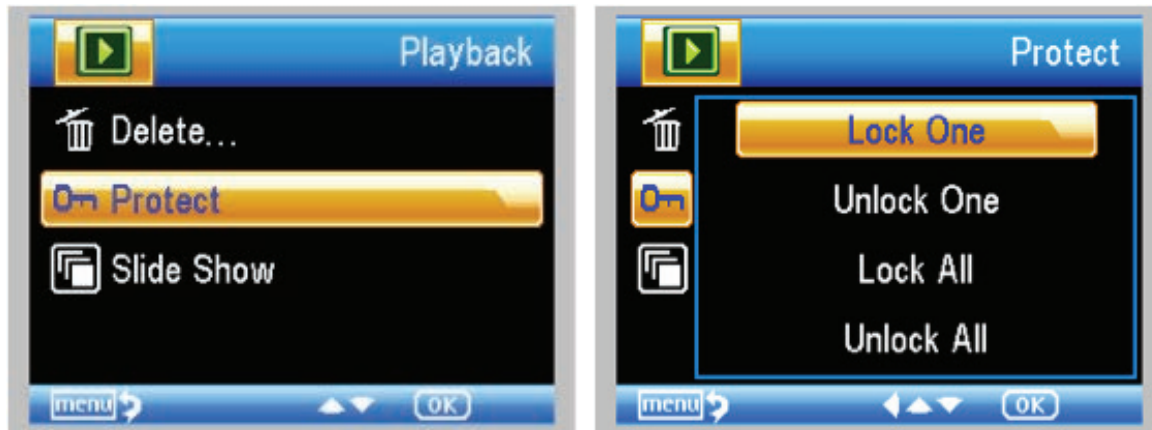

3. SLIDE SHOW:


This choice creates a slide show using all the image files on the MicroSD card. Select Slide Show, and then choose the interval between each image.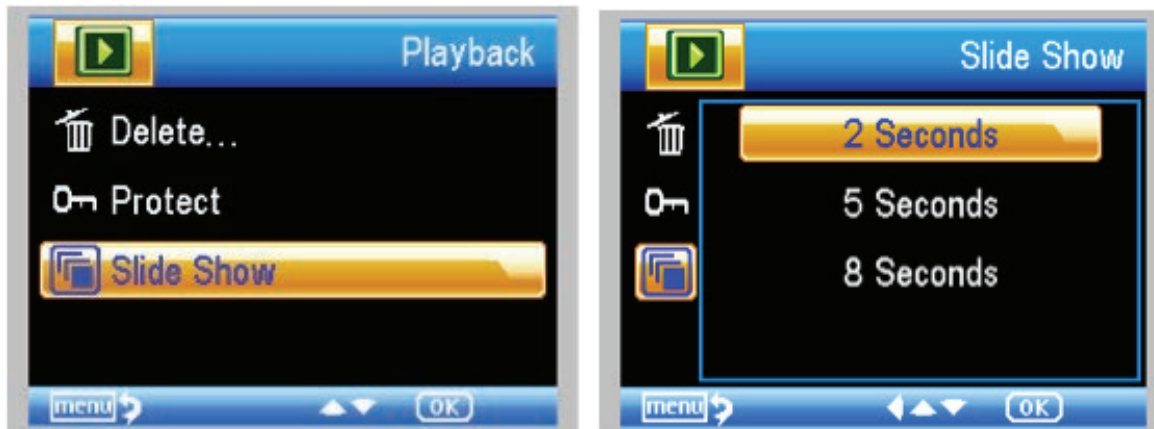

FILTERS:


When on the main screen or in viewing mode, press the MODE/UP button **M** ▲ to scroll between:

- A. Color Image
- B. Negative Image
- C. Black and White Image

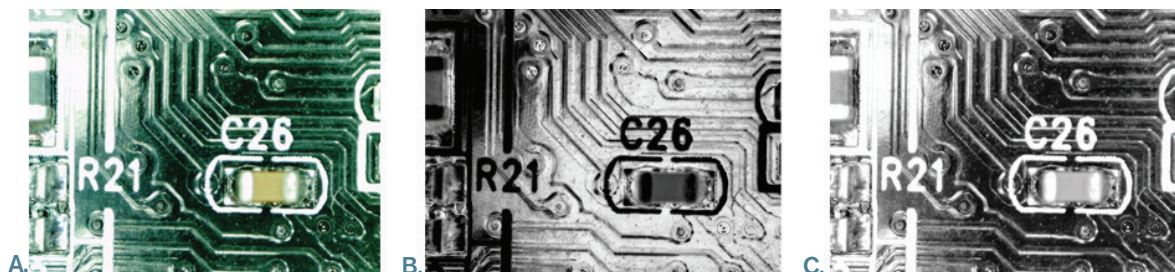DIGITAL ZOOM:

- + Press the Zoom Button up to zoom in, and down to zoom out.
- + The Zoom has a range from 1.0x to 4.0x

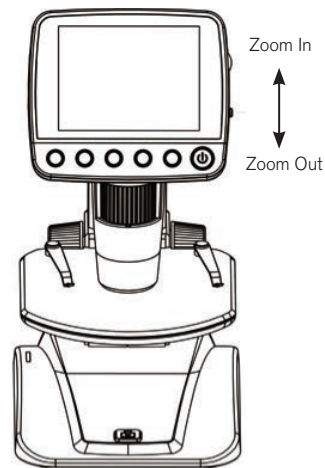6. SOFTWARE MENU INTRODUCTIONS

MAIN MENU 1 :

- + Press  SETTINGS button to enter Main Menu 1.
- + The menu choices will vary depending on whether you are in still image or video mode.
- + When in still image mode, you will see:
 1. Resolution
 2. Time Lapse
 3. Quality
 4. Sharpness

- + There is only one setting in video mode:

1. Resolution

RESOLUTION SETTING:

- + In still image mode, choices are:
 - 12.0 MP (interpolated*)
 - 9.0 MP (interpolated*)
 - 5.0 MP
 - 3.0 MP
 - 1.3 MP
 - VGA

+ In video mode, choices are:

- VGA
- QVGA

TIME LAPSE SETTING

+ FRAMES: In this window, you will set the number of frames you want to take.

NOTE: The number above represents 32 frames.

+ INTERVAL: In this window the interval between shots is set.

- Once you have made your selections, press OK to confirm and press the Menu/Setup button to go back to preview mode.
- When you press the Capture Button to begin your time lapse, you will notice a new icon  is displayed in the upper left corner of the screen.
- When all the frames are taken, the InfiniView will go back to the preview screen.
- To stop the auto capture before it completes, press the Capture Button once.

-NOTE: The TIME LAPSE setting only works in Still Image mode.

NOTE: The number above represents 15 seconds.

QUALITY

The Quality settings determine the level of compression for the image you are about to take. The more compression, the smaller the file size, but the less detail in the image.

+ There are 3 choices in the Quality setting window:

1. Super Fine (largest files, least compressed)
2. Fine
3. Normal (smallest files, most compressed)

SHARPNESS

Sharpness increases the differences the imaging process detects between the edges of the pixels. There are no rules when it comes to the sharpness settings. Experiment to find out which works best for your specimens.

+ The Sharpness choices are:

1. Hard
2. Normal
3. Soft

MAIN MENU 2

While in MAIN MENU 1, press the left button to enter MAIN MENU 2. The choices here are:

1. Language
2. Power Save
3. Frequency
4. Date Stamp
5. TV Out
6. Date/Time
7. Format SD
8. System Reset
9. Version

LANGUAGE

InfiniView may be used in 9 languages:

1. English
2. German
3. French
4. Spanish
5. Italian
6. Portuguese
7. Japanese
8. Simple Chinese
9. Traditional Chinese

+ Scroll to your preferred language and press OK to select it.

POWER SAVE

Power Save allows you to conserve battery power by setting the unit to turn off automatically. The choices for times are:

1. OFF: Power Save feature is off; unit will not turn off automatically
2. 5 min.: Unit will turn off after 5 minutes of non-use
3. 10 Min.: Unit will turn off after 10 minutes of non-use
4. 15 Min.: Unit will turn off after 15 minutes of non-use
5. 30 Min.: Unit will turn off after 30 minutes of non-use

FREQUENCY

This setting has to do with your TV's refresh rate. For USA (NTSC), use 60Hz; for others, use 50Hz.

DATE STAMP

If you wish, you can display a date stamp on image files (but not on video files).

+ The choices are:

1. DATE/TIME: Displays both DATE and TIME on the saved image file
2. DATE: Displays only the date on the image file
3. OFF: Will NOT display any DATE/TIME information


NOTE: The DATE and/or TIME will NOT be displayed in the preview mode—only on the recorded image file.

TV OUT

Set your TV out setting to either NTSC or PAL, depending on your location.

1. Choose NTSC for:
 - North America
 - South America (excluding Brazil, Argentina, Uruguay and French Guiana)
 - Myanmar, South Korea, Taiwan, Japan and the Philippines
2. Choose PAL for:
 - All countries NOT listed above

NOTE: See map below.

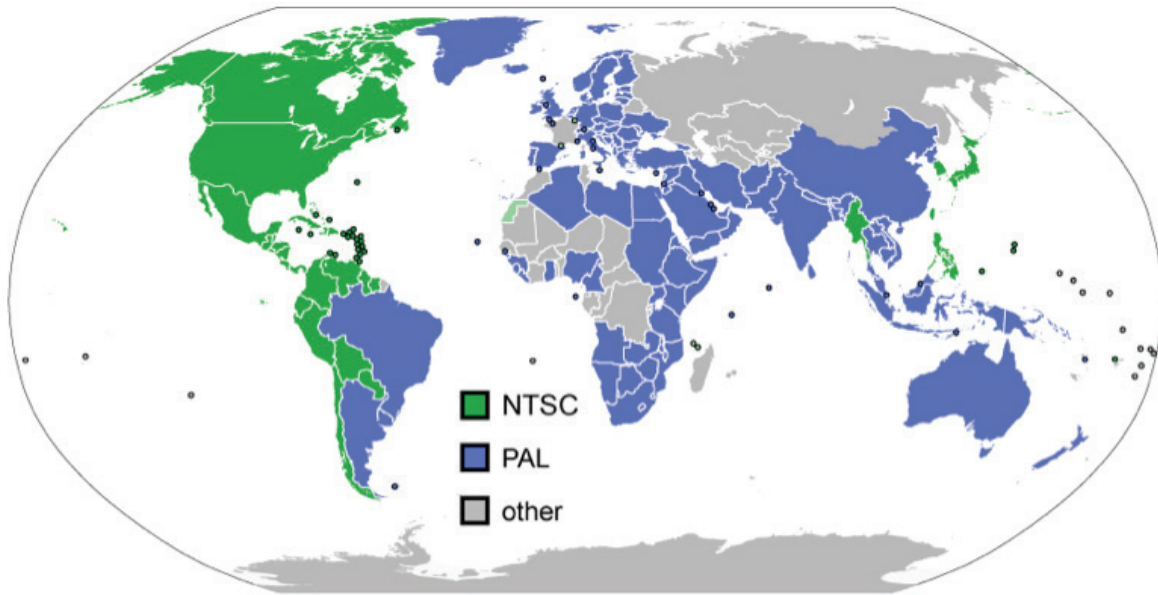DATE/TIME

- + Here you can set the correct Date and Time
- + You can also choose from the following date display options:
 - YY/MM/DD
 - DD/MM/YY
 - MM/DD/YY


SD CARD

From this screen, you can format your MicroSD card. You should format your MicroSD before using it for the first time.

NOTE: During formatting, ALL files will be erased from the card.

SYSTEM RESET

This screen allows you to reset ALL the settings to the factory default settings.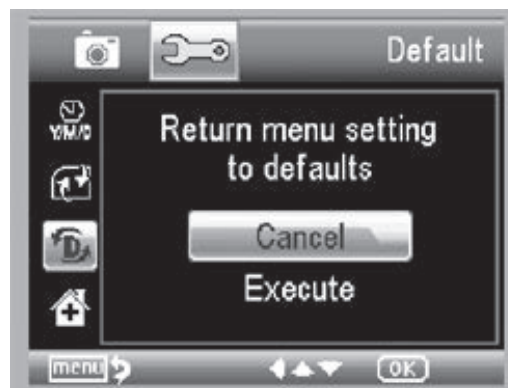


VERSION

The Version screen will display information on your firmware version. You cannot actually change any setting here.


TRANSFER FILES TO COMPUTER

- + Connect your InfiniView to your computer via USB 2.0
- + On the LCD screen, choose “Upload Images”, using the OK button.
- + Your computer will copy the files on your MicroSD card to your hard drive.
- + When the transfer is complete, you can open the files on your computer.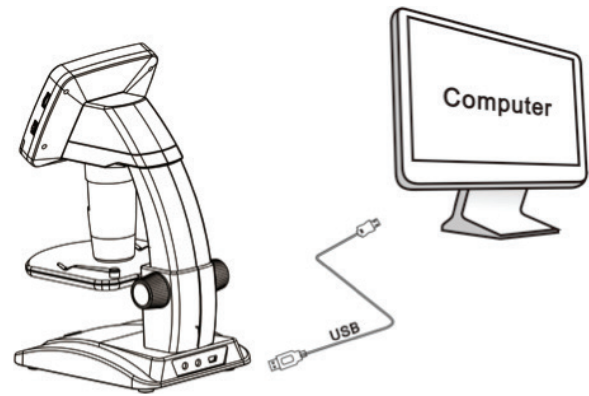

CONNECT TO TV

- + Insert the 1/8-inch plug on the end of the supplied video cable into the InfiniView AV OUT port located on the back of the base.
 - + Insert the yellow RCA plug on the video cable into one of your television's yellow video input connectors.
- NOTE:** Refer to your television's manual to locate your TV's video connector.
- + Turn on the television. (Be sure your television is tuned to the appropriate input channel.)
 - + You should now see an image of the specimen you are viewing on your television screen.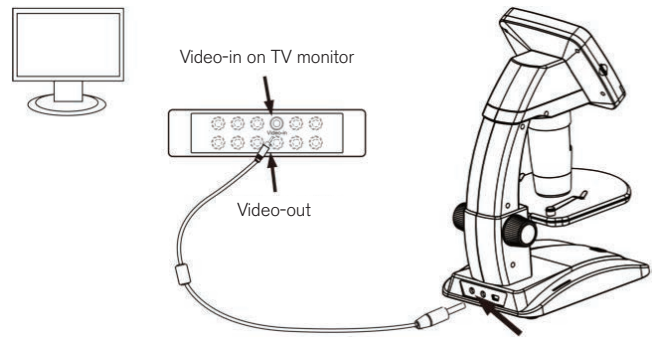

USING THE CELESTRON PORTABLE CAPTURE PRO SOFTWARE

- + Download and install the latest Portable Capture Pro software from the product page at celestron.com.
- + Connect the InfiniView to your computer as described above.
- + Choose “Preview on Computer” Mode from the 2 choices on the LCD screen.
- + Open the Celestron Portable CapturePro Software on your PC.

NOTE: Once the software is launched, the LCD screen will go blank and you will begin controlling the microscope with the computer. However, you will still use the InfiniView to control the features below.

NOTE: When you connect to a computer via USB, the lithium battery will charge.

7. USING YOUR MICROSCOPE

A. CONNECTING

+ Connect the microscope to your computer via the USB 2.0 cable.


+ Click on the Micro CapturePro desktop icon to launch the software.

+ Your microscope should automatically connect and begin streaming images.

NOTE: If the microscope is not connected to the USB port, the following error message will be displayed: "No device detected. Please connect your microscope directly to an open USB port."

1. Illuminator Brightness
2. Focus Wheel
3. Stage Adjustment Wheel
4. Stage Lock Wheel
5. Power button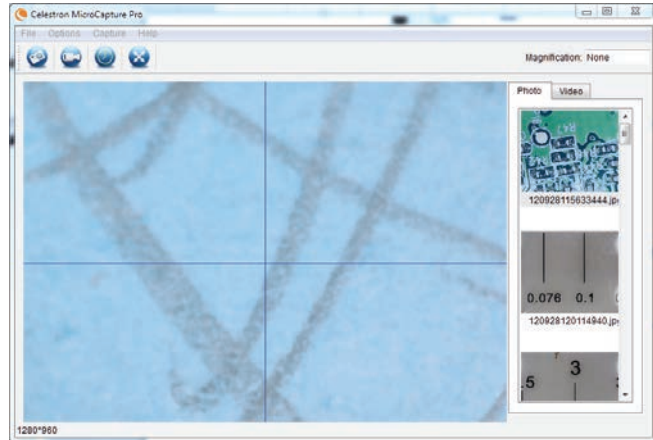

8. USING THE SOFTWARE

A. CAPTURE A PHOTO/VIDEO

+ To capture a photo, click on the Capture menu and scroll down to Photo.

+ To begin recording video, click on the Capture menu and scroll down to Video. Repeat to stop recording.

B. OTHER MENU ITEMS

FILE: Under the File menu item, you will find:

1. Photos Directory: Directs you to designate a folder on your hard drive for photo storage.
2. Videos Directory: Directs you to designate a folder on your hard drive for video storage.
3. Exit: Closes the Micro CapturePro software.

OPTIONS: Under the Options menu item, you will find:

1. Resolution: Sets the sensor image resolution.
Choose from 2 resolution settings:

- 320 x 240



- 640 x 480

NOTE: When connected to a PC, InfiniView cannot capture high-resolution images, but can be used for demonstration and streaming.
To take advantage of full 5 MP resolution, unplug InfiniView from the computer.


2. Date/Time: Turns the Date and Time stamp on or off.
3. Language: Changes the language setting. Choose from 12 languages:
 - English, German, French, Spanish, Russian, Italian, Portuguese, Dutch, Polish, Japanese, Korean and Chinese
4. X,Y Guide: Turns the reticle guide on the preview screen on or off.
5. Full Screen Mode: Enters full screen mode. Press ESC to exit.

ICONS: Along with the menu items, there are icons which mean the following:


- A. Take Snap Shot
- B. Start Video. Stop Video
- C. Set Time Lapse Intervals
- D. Full Screen Mode (press ESC to go back to main screen)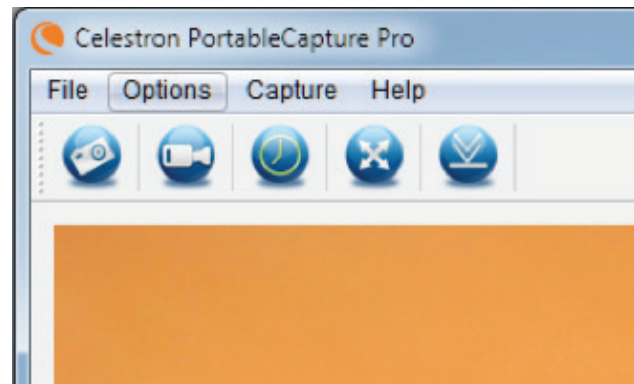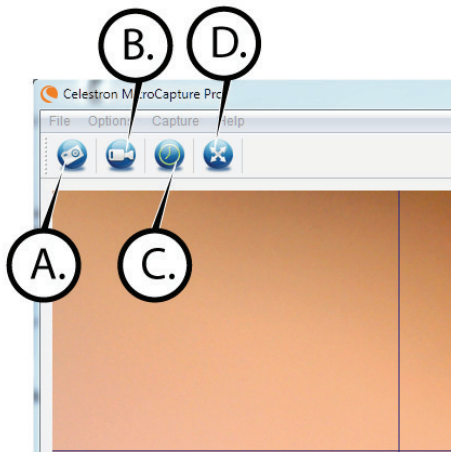


C. CALIBRATION AND MEASUREMENT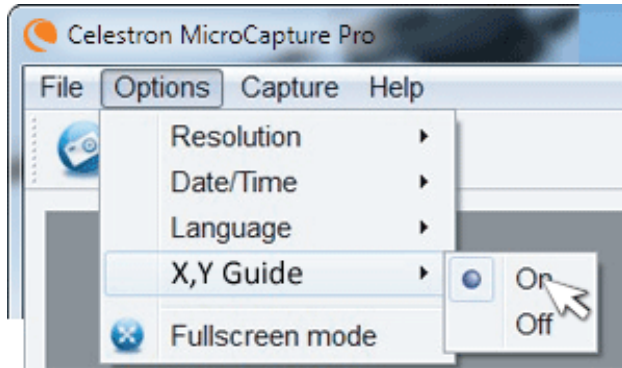


- + The Celestron Micro CapturePro software includes a measurement feature. To use this feature, you must first perform a measurement calibration.
- + Make sure the X,Y Guide is set to "On" by selecting it in the Options Menu.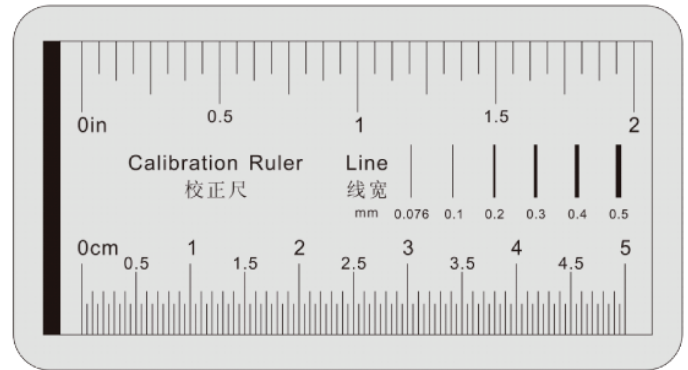

- + Set the height of the microscope and focus on your specimen. Then, replace your specimen with the calibration ruler.
- + Re-adjust the focus and align the calibration ruler so that the lines on the ruler are parallel with the vertical line on the X,Y Guide.

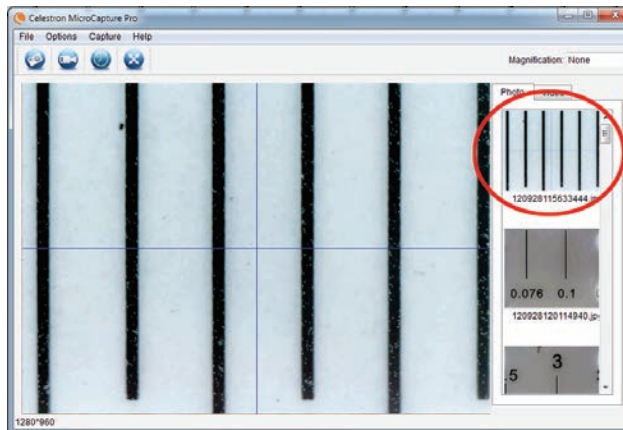
- + Take a still image. The image will be displayed in the column on the right side of the screen.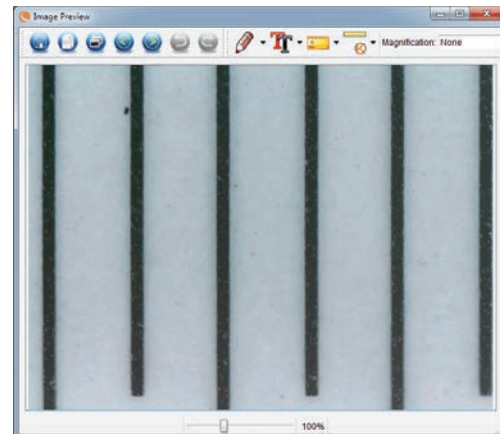

- + Double click the saved image to open it in the Image Preview screen.

- + Click on the Calibration icon 
- + Click and drag a line across a "known value" on the ruler image. Try to keep the line as perpendicular to the lines on the ruler as possible.
- + The "Adjust magnification" window will pop up as shown here.
- + In the box "Actual Dimension", enter the "known value" from the calibration ruler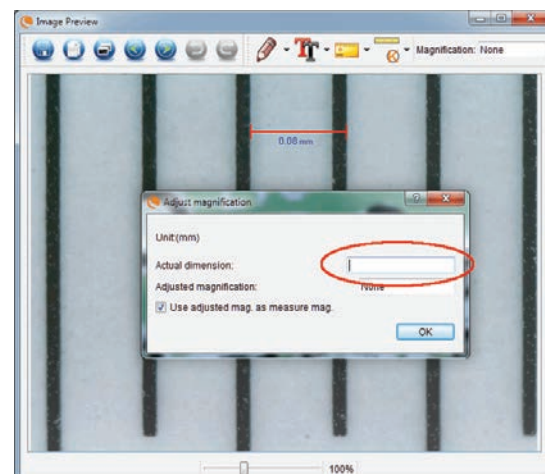

+ To ensure that calibration was successful, go back to the Image Capture screen.

+ Click the ruler icon  and from the dropdown menu select the option "Any Angle Line."

+ Click and drag a line across the same "known value" as before.

+ If the calibration was successful, the distance you measured should be exact to the "known value."

+ To capture images without changing the height or magnification, click on the Calibration icon. In the dropdown menu, select "Set Magnification," then "Set Capture Magnification."

+ A pop-up window will appear. Type in the magnification shown in the upper right corner of the main window.

D. MEASURING TOOLS




You will notice in the measuring drop down menu there are many tools. Here is the list of measurements you can take:

MAIN

1. Any Angle: Measures from point A to point B in a straight line.
2. Continuous Line: Measures the total distance between multiple points. Start by measuring a straight line from point A to point B. Then, click and drag to continue measuring from point B to point C. (Can be repeated as many times as you like.)
3. Radius Circle: Measures a circle's radius. Starting at the center, drag the cursor out to any edge.
4. Diameter Circle: Measures a circle's diameter. Starting at one outside edge, drag the cursor across to the opposite edge.
5. Three Point Angle: Measures an angle between two points, starting from a radius point or an intersection point. Click on point A, (radius point) drag and click on point B, then click on point C. The angle will be drawn and the angle degree will be displayed.

SETTINGS

1. Line Properties: Changes the settings for line type, width, size and color.
2. Text Properties: Changes the settings for font type, size, style and color.
3. Set Units: Changes the units setting. Choose from millimeters, centimeters, pixels or inches.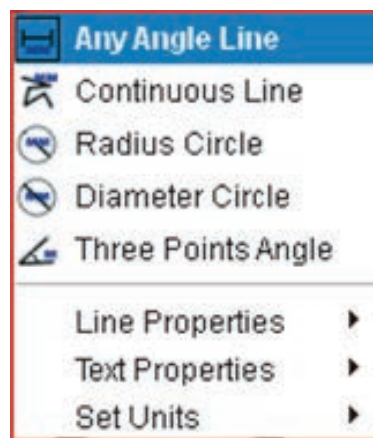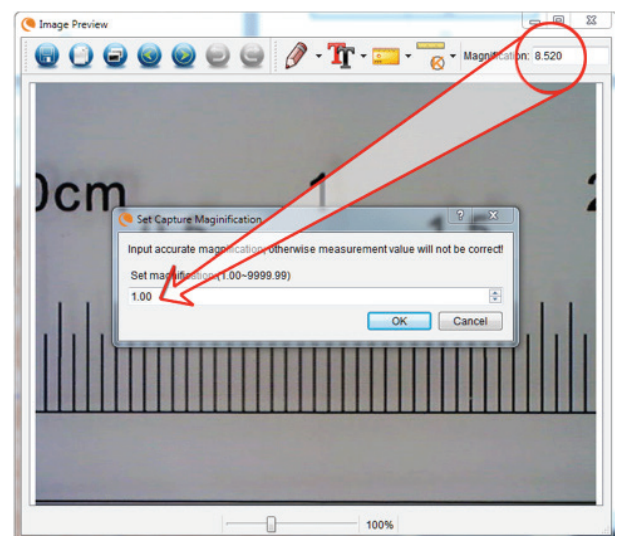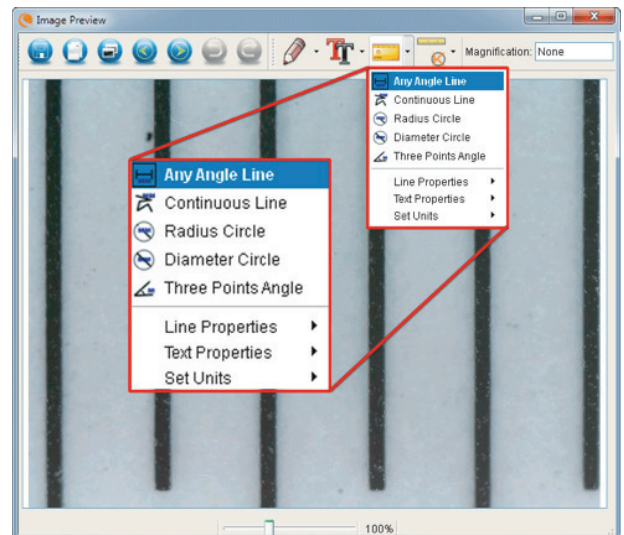

E. TAKING NOTES AND MARKING THE IMAGE

Micro CapturePro includes functionality to mark, draw on, or add notes to saved images. This can be handy for quick note taking and pointing out features on your specimen image.

1. MARKING


- + To use this feature, click the Draw icon  in the Image Preview window.
- + The complete choices in the Draw dropdown menu are:

MAIN

1. Line: Draws a straight line on the image.
2. Arrow: Draws a line with an arrow on the image.
3. Freehand: Allows user to draw freehand. Good for quick, handwritten notes.
4. Rectangle: Draws a box. Good for calling out features on the image or framing text.


SETTINGS

1. Line Properties: Allows for setting the line type, size and color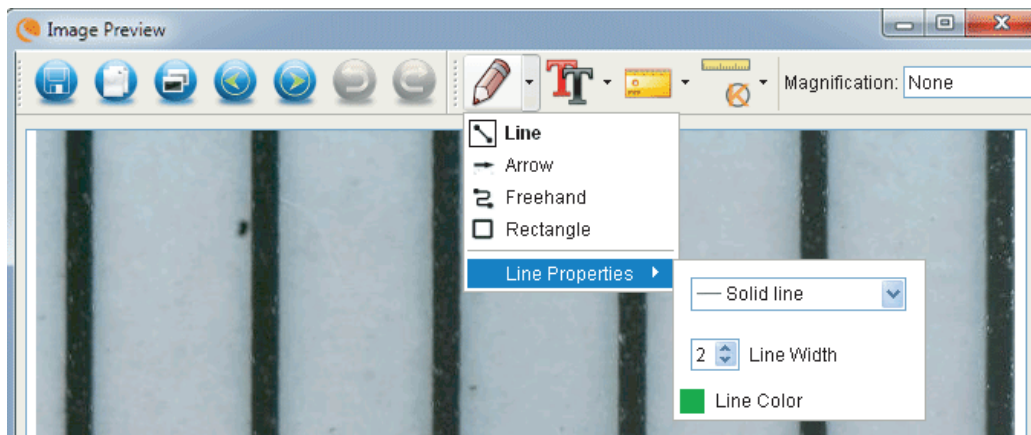


2. ADD TEXT

- + To type notes on the image, select the Type icon .
- + In the dropdown menu, you can choose the font type, size and color.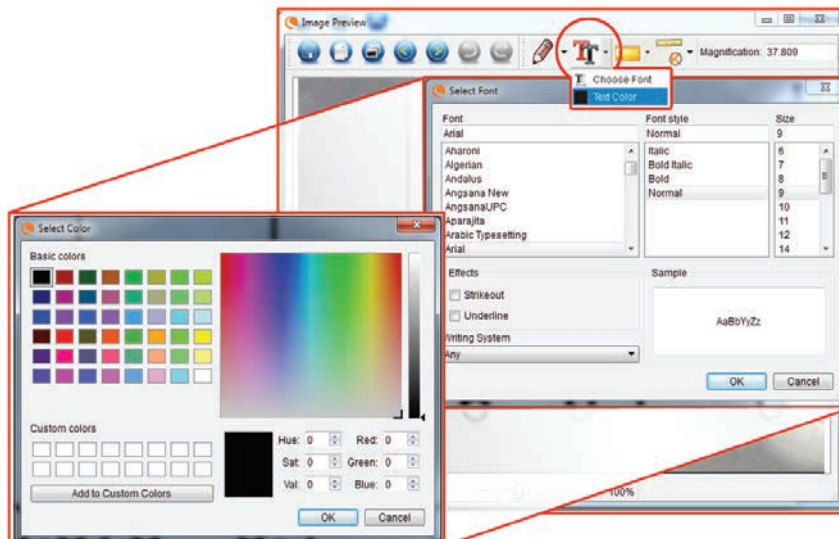

F. IMAGE PREVIEW ICONS EXPLAINED

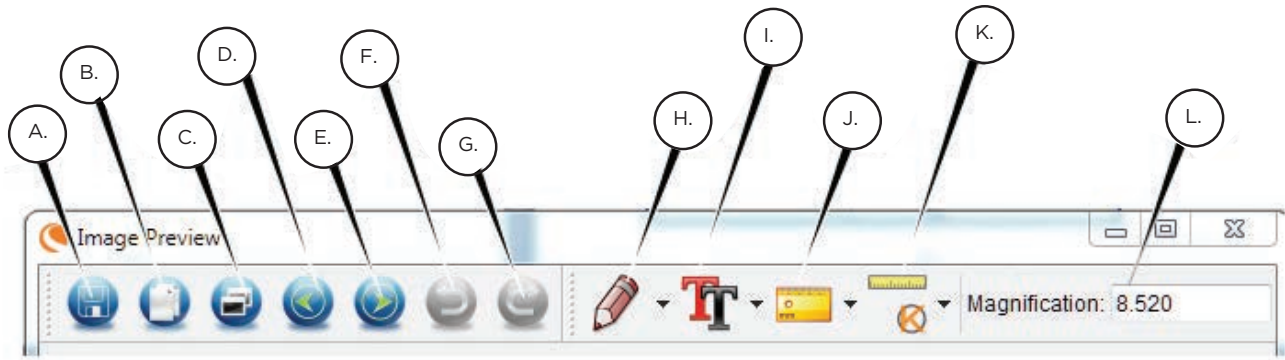- A. Save: Save your file
- B. Copy to clipboard: Copy to open in other program
- C. Copy: Save as a copy
- D. Previous: Takes you back to the previous image in save folder
- E. Next: Takes you forward to the next image in save folder
- F. Undo: Takes you back one step
- G. Redo: Takes you forward one step
- H. Draw
- I. Text/Notes
- J. Measure
- K. Calibration
- L. Current Magnification (based on the latest calibration)

9. MULTI-PLUG

If you have purchased your InfiniView outside of North America (model 44361), you will notice a multi-plug AC adapter. There are 3 interchangeable plug heads that will fit outlets in Europe, the United Kingdom, Hong Kong and China.

To change the plug head:

- Grasp the main plug housing with one hand and the plug head of choice with the other.
- Press the small button release on the plug housing (circled in below image).
- Locate and interlock the tabs on the plug head with the grooves on the plug housing.
- Rotate the plug head clockwise 90° until you hear a "click" as the plug head locks into place.

To remove the plug head, press the small button release and rotate the plug head counterclockwise until it releases.

10. CARE AND MAINTENANCE

Your microscope with digital camera is a precision optical instrument and should be treated with care at all times. Follow these care and maintenance suggestions and your microscope will need very little maintenance throughout its lifetime.

- + Store your microscope in a dry, clean place.
- + Never use your microscope where operation of such devices is restricted. Improper use creates the risk of serious accident.
- + Do not try to access the internal components of your microscope or attempt to modify it in any way. Only authorized technicians should perform maintenance or repairs.
- + Keep your microscope away from water and other liquids. Never use it in the rain or snow. Moisture creates the danger of fire and electric shock.
- + To clean dust or smudges from the LCD screen or other optical component, first blow the dust off with a camel's hair brush or air blower. Then use the included dry, microfiber cloth to wipe the surface.
- + If the exterior surfaces of your microscope need cleaning, wipe with a moist cloth.

11. TROUBLESHOOTING

PROBLEMS	SOLUTION
The unit does not turn on	Check the battery power. If necessary, recharge by plugging into AC power
The screen is blank	<ul style="list-style-type: none"> + Ensure that the power is on + Check the battery power and recharge the battery if necessary + Verify that the object you want to magnify is positioned correctly under the lens
There are smudges or blemishes on the screen	Clean the screen and lens carefully with the included microfiber cloth
Microscope is connected to a television, but no image appears on the television screen	<ul style="list-style-type: none"> + Ensure that the television is on. + Switch to the appropriate video input channel on your television (this will vary by manufacturer) + Ensure that the video cable is securely connected to the microscope's Video Out port and plugged into the correct video input on your television
The battery indicator LED is not illuminating, even though the power adapter is connected and plugged into a wall socket -or- The battery is not charging	<p>Make sure the cable is plugged into the proper port.</p> <p>Check that the outlet has power and is not malfunctioning. If the outlet is connected to a wall switch, make sure that the switch is on</p>

12. WARRANTY

Your InfiniView LCD Microscope has a two year limited warranty. Please visit the Celestron website for detailed information on all Celestron microscopes: celestron.com

© 2020 Celestron • All rights reserved.
celestron.com/pages/technical-support
 2835 Columbia Street • Torrance, CA 90503 USA

FCC Note: This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- + Reorient or relocate the receiving antenna.
- + Increase the separation between the equipment and receiver.
- + Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- + Consult the dealer or an experienced radio/TV technician for help.

This product is designed and intended for use by those 14 years of age and older. Product design and specifications are subject to change without prior notification.

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