HD DIGITAL **MICROSCOPE IMAGER**

QUICK SET-UP GUIDE

MODEL #44422

INTRODUCTION

Congratulations on your purchase of the Celestron HD Digital Microscope Imager. Your Imager is a precision optical instrument made of high quality materials to ensure durability and long life, and designed to give you a lifetime of use with minimal maintenance. It is miniature in size – 2.5" (64mm) long x 0.9" (22mm) in diameter and weighs just 1.5oz (42g) but it is powerful in its application with a 5MP sensor digital CMOS sensor.

The HD Digital Microscope Imager will work with virtually any microscope (biological and stereo) as long as the eyepiece barrels are universal sizes 23mm and 30mm or similar. Basically, you replace the eyepiece in the microscope with the Imager.

Before attempting to use your Imager, please read through the instructions to familiarize yourself with the functions and operations to make usage easier for you.

Compatibility with Windows and Mac operating systems allows for high resolution images and video to be taken, stored on your computer, saved and shared.

NOTE: The Mac software only supports a maximum 2MP resolution.

USING YOU HD DIGITAL MICROSCOPE IMAGER

1. Remove the HD Imager and the collar from the box.
2. Attach the adjustment collar to the HD Imager using the set screws to secure.
3. Remove the eyepiece from your microscope and replace it with the assemble imager and collar.
4. Install the software from the product page at celestron.com, on to your Windows or Mac PC.
NOTE: Mac operating systems are supported, but can only achieve a maximum of 2MP resolution when taking still images.
5. Once the software is installed, connect your imager by plugging the attached USB cable into an open USB port on your computer.
6. Once the driver is recognized by your system, the streaming image should appear in your software window.
7. Focus the microscope as you would normally.

APPROXIMATING THE FINAL MAGNIFICATION

The approximate power is determined by the Imager power multiplied by the objective lens power. The Imager has a 30x magnifying lens built in. As an example, if the objective lens being used is 4x, then the power you will be capturing for snapshots or video will be 120x. If using a 40x objective, then the power will be 1200x, etc.

ROTATING THE IMAGER

When you are looking at the image on your computer screen, you can rotate the position by rotating the Imager to the position you prefer.

USING THE CELESTRON IMAGER HD SOFTWARE

The software is very easy to use and should be very intuitive. there are many features that you can control however, so if you wish to delve deeper into the Celestron Imager HD software, click the Manual icon in the software  or download the software manual from the product page at celestron.com.

YOU ARE NOW READY TO BEGIN YOUR DIGITAL ADVENTURE!

CARE, MAINTENANCE, AND WARRANTY

Your Imager is a precision optical instrument and should be treated with care at all times. Follow these care and maintenance suggestions and your Imager will need very little maintenance throughout its lifetime.

- Store the Imager in a dry and clean place.
- Be very careful if using your imager and microscope in direct sun light to prevent damage to the Imager or your eyes.
- Never point the Imager towards the sun or the sensor can be damaged and cease working.
- To clean the housing, use a damp cloth and dry immediately.
- Use a good lens tissue to clean the clear glass end (optical surface) of the Imager.
- If needed, blow off dust with a camel's hair brush or air blower off the optical surfaces.
- To clean fingerprints off of optical surfaces, use a lens cleaning agent and lens tissue available at most photo outlets and when cleaning do not rub in circles as this may cause streaks or scratches to occur.
- Never disassemble or clean internal optical surfaces. This should be done by qualified technicians at the factory or other authorized repair facilities.
- Do not disassemble or take apart the Imager or damage can occur.

WARRANTY

Your Imager has a two year limited warranty. Please see the Celestron website for detailed information at www.celestron.com.

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FCC Note: This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- + Reorient or relocate the receiving antenna.
- + Increase the separation between the equipment and receiver.
- + Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- + Consult the dealer or an experienced radio/TV technician for help.

©2020 Celestron • All rights reserved.

celestron.com/pages/technical-support

2835 Columbia Street • Torrance, CA 90503 USA

This product is designed and intended for use by those 14 years of age and older. Product design and specifications are subject to change without prior notification. Version 2.

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